



Please note that booster heaters are available for all models except P1-C25P and DC30-K20 which will reduce heating times by as much as 60%

Reactor	Time to Cool to -60	Time to Cool to -20	Time to Heat to +100	Time to Heat to +90
1 Liter	P2-CT80L: 1 hour and 40 minutes	DC30-K20: 55 minutes P1-C25P: 55 minutes P1-C35P: 70 minutes P1-C50P: 35 minutes	DC30-K20: 25 minutes P1-C25P: 25 minutes P1-C35P: 40 minutes P1-C50P: 25 minutes	P2-CT80L: 1 hour
2 Liter	P2-CT80L: 1 hour and 50 minutes	DC30-K20: 1 hour and 10 minutes P1-C25P: 1 hour and 10 minutes P1-C35P: 1 hour and 25 minutes P1-C50P: 40 minutes	DC30-K20: 30 minutes P1-C25P: 30 minutes P1-C35P: 45 minutes P1-C50P: 28 minutes	P2-CT80L: 65 minutes
3 Liter	P2-CT80L: 2 hours	DC30-K20: 1 hour and 23 minutes P1-C25P: 1 hour and 23 minutes P1-C35P: 1 hour and 35 minutes P1-C50P: 45 minutes	DC30-K20: 37 minutes P1-C25P: 37 minutes P1-C35P: 52 minutes P1-C50P: 31 minutes	P2-CT80L: 1 hour and 15 minutes
4 Liter	P2-CT90L: 1 hour and 15 minutes	DC30-K20: 1 hour and 40 minutes P1-C25P: 1 hour and 40 minutes P1-C35P: 1 hour and 50 minutes P1-C50P: 50 minutes	DC30-K20: 45 minutes P1-C25P: 45 minutes P1-C35P: 1 hour P1-C50P: 35 minutes	P2-CT90L: 47 minutes
5 liter	P2-CT90L: 1 hour and 20 minutes P2-CT90W: 1 hour and 10 minutes	P1-C35P: 2 hours P1-C50P: 55 minutes P1-C30P: 1 hour and 15 minutes	P1-C35P: 65 minutes P1-C50P: 39 minutes P1-C30P: 50 minutes	P2-CT90L: 50 minutes P2-CT90W: 50 minutes
10 Liter	P2-CT90L: 1 hour and 45 minutes P2-CT90W: 1 hour and 30 minutes	P1-C50P: 1 hour and 20 minutes P1-C30P: 1 hour and 45 minutes P2-CT50L: 45 minutes P2-CT50W: 23 minutes	P1-C50P: 1 hour P1-C30P: 1 hour and 10 minutes	P2-CT50L: 1 hour and 25 minutes P2-CT50W: 58 minutes P2-CT90L: 68 minutes P2-CT90W: 50 minutes
15 Liter	P2-CT90L: 2 hour and 10 minutes P2-CT90W: 1 hour and 55 minutes	P1-C50P: 1 hour and 45 minutes P1-C30P: 2 hours and 10 minutes P2-CT50L: 55 minutes P2-CT50W: 26 minutes	P1-C50P: 1 hour and 18 minutes P1-C30P: 1 hour and 28 minutes	P2-CT50L: 1 hour and 40 minutes P2-CT50W: 1 hour and 10 minutes P2-CT90L: 1 hour and 22 minutes P2-CT90L: 1 hour and 22 minutes
20 Liter	P2-CT50L: 1 hour and 5 minutes P2-CT50W: 32 minutes P2-CT90L: 2 hours and 30 minutes P2-CT90W: 2 hours and 15 minutes	P1-C50P: 2 hours and 5 minutes	P1-C50P: 1 hour and 35 minutes P2-CT50L: 1 hour and 55 minutes P2-CT50W: 1 hour and 18 minutes	P2-CT90L: 1 hour and 35 minutes P2-CT90W: 1 hour and 35 minutes
30 Liter	P2-CT50L: 1 hour and 25 minutes P2-CT50W: 42 minutes P2-CT90W: 3 hours		P2-CT50L: 2 hours and 35 minutes P2-CT50W: 1 hour and 40 minutes	P2-CT90W: 2 hours and 5 minutes
50 Liter	P2-CT50W: 1 hour		P2-CT50W: 2 hours and 20 minutes	